

Firmware Updating Steps:

1. Unzip the downloaded folder (you may need to close this note first).
2. Open the "VCPLUS" Application.
3. Plug in the VCPlus 8 Valve device to power.
4. Turn on the power switch for the VCPlus 8 Valve device.
5. Plug in the USB Type C cable into the VCPlus 8 Valve device and into your computer.

(You can use either USB Type-C to Type-A or Type-C to Type-C to connect the VCPlus device to the computer).
6. Press Settings>Update Firmware>Load File.
7. Search for the unzipped downloaded "VC8_5468_8C_v2.3" file and double click it.
8. Press the "Program" button and wait for the "Firmware Upgraded Sucessfully!" pop up.
9. You can now close the Firmware updating window(s).